



Mega-Pure® Accessories ACS



Automatic Collection System (ACS)

Automatic Collection = No Hassle!

- All glass design with ACS
- ACS designed to collect water from Mega-Pure stills
- 45 liter capacity

Product Description

- The MEGA-PURE Automatic Collection System is designed to collect water from the MEGA-PURE stills and control their operation.
- 45 liter capacity.
- Wall or bench mounted.
- All glass design.
- Glass inlet and outlet tubing.

Operation

- ACS will work with MEGA-PURE stills to store up to 45 liters of Type II distilled water.

- The ACS is preset to shut off the power to the heaters and the water supply to the MEGA-PURE still after the bottle has collected approximately 45 liters.
- When the purified water supply has been depleted to approximately 38 liters, the system will automatically restart the water still and refill the collection system.
- When setting up a MEGA-PURE still with the ACS on a bench, a 6-12" stand (depending on still model) is required. (This will ensure proper gravity feed of distilled water.)

PRODUCT SPECIFICATIONS

Model #	W	Dimensions Inches (cm)		Operating Weight Lb. (kg)	Shipping Weight Lb. (kg)
		H	D		
B440704	16.25 (41.3)	38 (96.5)	16.25 (41.3)	110 (49.9)	66 (29.9)

ORDERING INFORMATION

Model #	Power	Capacity liter	US List Price
B440704	None	45	\$1,922.00

ACCESSORIES

Model #	Where Used	US List Price
ACS		
440138	Flexible Tubing Adapter Kit —When placing ACS in not recommended set-up ¹	\$102.00
400634	Wall Mounting Bracket for ACS	\$227.00
440241	Adapter for ACS Washer Hook-up for Glassware Washers	\$212.00
D6832	ACS to NANOpure Infinity/Diamond Adapter	\$353.00

¹ Requires customer supplied 1/2" ID Teflon tubing.